

Analysis of Limestone for Flue Gas Desulfurization in a Power Plant

ARL OPTIM'X Simultaneous-Sequential Wavelength Dispersive X-Ray Fluorescence Spectrometer

Key Words

- ARL OPTIM'X
- Lime/limestone
- Power plant
- WDXRF
- X-Ray Fluorescence


Industry background

What is flue gas desulfurization?

A flue gas desulfurization (FGD) processing unit, commonly referred to as a scrubber, removes sulfur dioxide (SO₂) from the exhaust flue gases in power plants that burn coal; airborne SO₂ forms acid rain. As environmental emissions regulations for sulfur dioxide tighten in countries worldwide, technologies for SO₂ removal and associated analysis continue to advance.

How does a scrubber work?

During the combustion process the sulfur in the coal combines with oxygen in the air to form SO₂. To remove the SO₂, the flue exhaust from a coal-fired power plant is commonly bubbled through a mixture of lime or limestone and water. The resultant reaction typically captures 95 % or more of the SO₂. Approximately 80-85 % of worldwide FGD units installed in power plants use wet limestone scrubbing vs. other technologies.

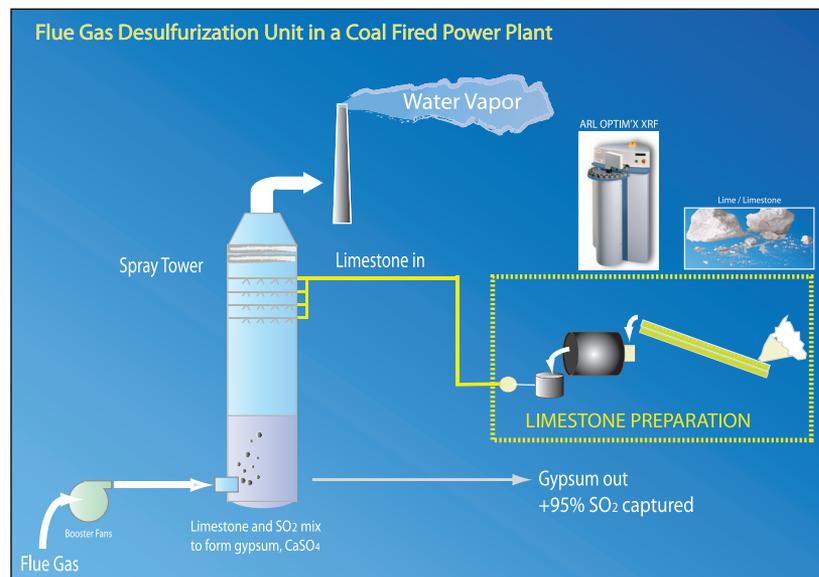

What happens to the SO₂ captured in a scrubber?

The captured SO₂ combines with the lime or limestone slurry (Ca(OH)₂ or CaCO₃) to form several byproducts, primarily calcium sulfate (CaSO₄) commonly called gypsum. Gypsum is recyclable or marketable for many industrial uses; the gypsum created by a power plant that is not reused is disposed of in permitted landfills along with scrubber byproducts that are not reusable.

XRF application

The acceptability of the lime or limestone used for scrubbing may be determined through element/oxide analysis, a perfect application for XRF. Typical ranges and reproducibility of interest for limestone would be:

Element/Oxide	Concentration	Required Std. Dev
CaCO ₃	> 95.0 %	0.10
SiO ₂	< 3.0 %	0.05
MgCO ₃	< 2.0 %	0.05
Al ₂ O ₃	< 1.5 %	0.05
Fe ₂ O ₃	< 3.0 %	0.03

Instrumentation

A Thermo Scientific ARL OPTIM'X XRF spectrometer has been used to derive limits of detection and precision for the analysis of lime/limestone. The ARL OPTIM'X is a WDXRF system which provides superior resolution and light elements capability. It is fitted with an air-cooled Rh end-window tube with thin Be window (0.075 mm) and has a maximum power of 50 Watts. Thanks to close coupling between the X-ray tube anode and the sample, the performance of the ARL OPTIM'X is equivalent to a 200 W conventional WDXRF instrument. The instrument can be equipped with the unique SmartGonio™, a series of Multichromators™ or both for various elements.



Instrument configuration

- ARL OPTIM'X analyzer – vacuum chamber, SmartGonio™, element range F to U
- 12-position sample changer
- OXSAS XRF analytical software

With an analysis time of only 10-40 seconds per element of interest, ARL OPTIM'X can easily satisfy the required repeatability with standard deviations as low as 0.001 to 0.03 depending on the element. Longer analysis times provide further improved results.

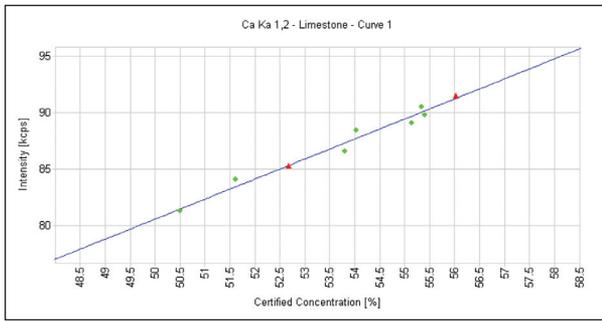


Figure 1: Calibration curve for calcium in limestone

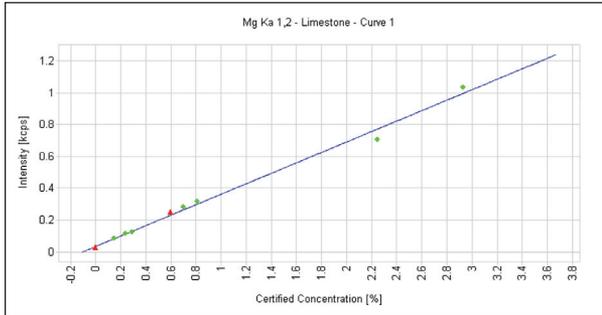


Figure 2: Calibration curve for magnesium in limestone

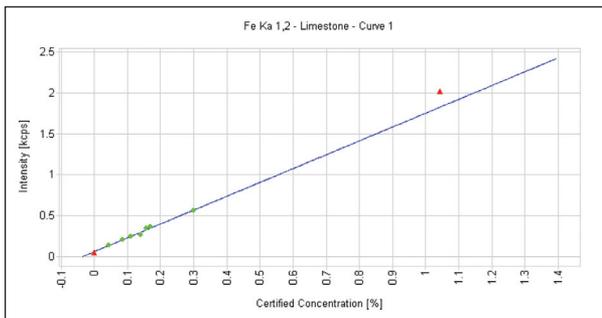


Figure 3: Calibration curve for iron in limestone

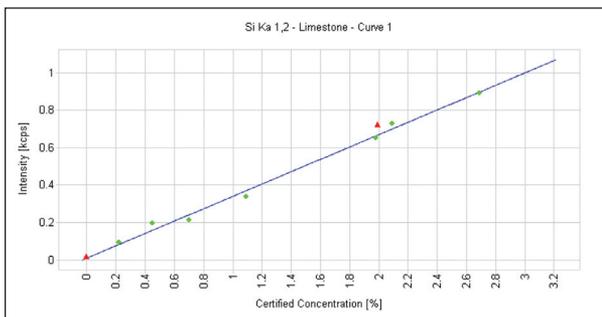


Figure 4: Calibration curve for silicon in limestone

Calibration and repeatability

A standard calibration curve has been developed on the ARL OPTIM'X for the five elements (Ca, Si, Mg, Al, Fe) in limestone as a pressed powder. The resulting calibrations for calcium, magnesium and iron are illustrated in Figures 1, 2 and 3.

A repeatability test was carried out by running a sample for ten repeat analyses with a counting time of two minutes per element. The results are summarized in Table 1.

Run #	CaCO ₃ (%)	SiO ₂ (%)	MgCO ₃ (%)	Al ₂ O ₃ (%)	Fe ₂ O ₃ (%)
Cell 1	95.56	2.10	1.60	0.49	0.26
Cell 2	95.35	2.11	1.60	0.49	0.26
Cell 3	95.59	2.10	1.60	0.50	0.26
Cell 4	95.67	2.09	1.61	0.50	0.26
Cell 5	95.31	2.12	1.58	0.50	0.26
Cell 6	95.43	2.10	1.60	0.50	0.26
Cell 7	95.57	2.12	1.59	0.50	0.26
Cell 8	95.41	2.10	1.61	0.49	0.26
Cell 9	95.48	2.08	1.59	0.50	0.26
Cell 10	95.47	2.10	1.59	0.50	0.26
Average	95.49	2.10	1.60	0.50	0.26
SD	0.114	0.012	0.011	0.005	0.002

Table 1: Repeatability for element oxides in pressed limestone powder

Conclusion

Excellent calibration curves and repeatability can be achieved with the ARL OPTIM'X to determine the suitability of limestone minerals for use in power plant scrubbers. The technique is both rapid and suitable for multiple oxides of interest at both high levels (approaching 100 %) and quite low levels (less than 0.25 %) of inclusion.

For more information on our X-ray product range and update on our latest developments, visit www.thermo.com/xray.

In addition to these offices, Thermo Fisher Scientific maintains a network of representative organizations throughout the world.

Africa-Other
+27 11 570 1840

Australia
+61 2 8844 9500

Austria
+43 1 333 50 34 0

Belgium
+32 2 482 30 30

Canada
+1 800 530 8447

China
+86 10 8419 3588

Denmark
+45 70 23 62 60

Europe-Other
+43 1 333 50 34 0

Finland/Norway/Sweden
+46 8 556 468 00

France
+33 1 60 92 48 00

Germany
+49 6103 408 1014

India
+91 22 6742 9434

Italy
+39 02 950 591

Japan
+81 45 453 9100

Latin America
+1 608 276 5659

Middle East
+43 1 333 50 34 0

Netherlands
+31 76 579 55 55

South Africa
+27 11 570 1840

Spain
+34 914 845 965

Switzerland
+41 21 694 71 11

UK
+44 1442 233555

USA
+1 800 532 4752

www.thermo.com



Thermo Fisher Scientific (Eublens) SARL, Switzerland is ISO certified.

AN41711_E 04/09C